



4. In the upper left corner of the LabVIEW screen,

click the Start button  which will change the button color to black .

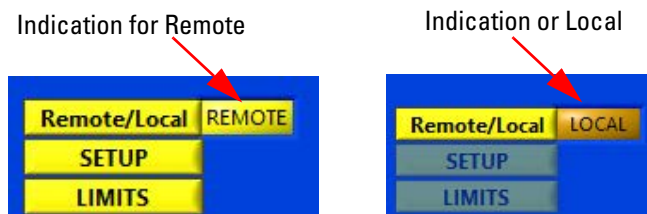


Figure 6: Remote/Local Indication

5. The box to the right of Remote/Local button should read either Remote or Local (See figure 6). Click on the Remote/Local button to toggle from Local to Remote for remote control. The controller screen should also show Remote on the lower left hand corner (See Figure 7).

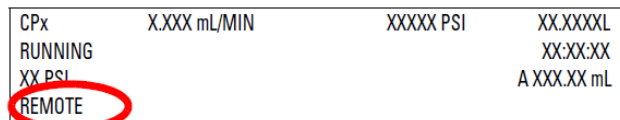


Figure 7: Controller display

Basic Settings and Commands

Some basic selections available in the sample program shown in Figure 8 (See Technical Bulletin TB39 for complete operation).

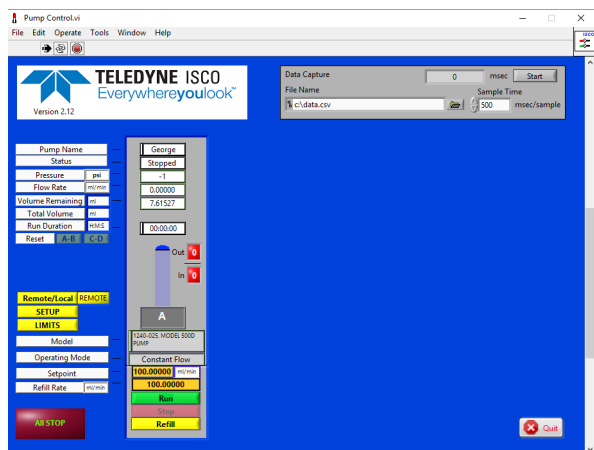


Figure 8: Example Program main screen

Pump Mode

To set Pump Mode, click on the SETUP button then select the MODE (Constant Flow, Constant Pressure, Continuous Const. Flow or Continuous Const. Pressure) for each pump under the MODE and PUMP designation (Figure 9).



Figure 9: Select operating mode

Set Points

On the main screen, click on the numbers to the right of the SETPOINT bar and enter the desired numeric value (See figure 10).

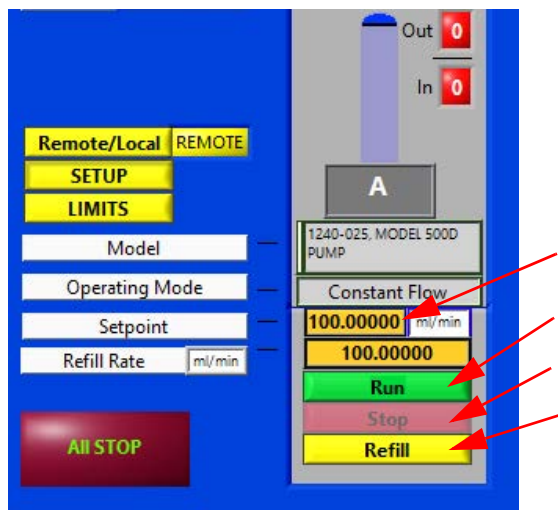


Figure 10: Set Points- Run, Stop, or Refill

Start/Stop

To Start, Refill or Stop each pump click either the RUN, STOP or REFILL buttons (See Figure 8).

Last modified May 12, 2023

Teledyne ISCO

P.O. Box 82531, Lincoln, Nebraska, 68501 USA
Toll-free: (800) 775-2965 • Phone: (402) 464-0231 • Fax: (402) 465-3001
E-mail: IscoService@teledyne.com

Teledyne ISCO is continually improving its products and reserves the right to change product specifications, replacement parts, schematics, and instructions without notice.



Use and Disclosure of Data

Information contained herein is classified as EAR99 under the U.S. Export Administration Regulations. Export, reexport or diversion contrary to U.S. law is prohibited.